

Pipette MethodCard

Operating Instructions



Pipette Performance Check
MethodCard

METTLER TOLEDO

1. Download EasyDirect - Pipette Asset Management (PAM) - Software



www.mt.com/SmartStand



Download & Install EasyDirect PAM

2. Connect RFID reader/writer to PC (see EasyDirect Manual)



PC with EasyDirect software



USB – RS Cable: 30091827



RFID Reader: 30215407

3. Define Method first, then write Method on Card (see EasyDirect Manual)



4. Hold card at XPE balance with RFID Reader (see XPE Balance Manual)



Follow instructions on the balance display to check the pipette

5. Edit or Delete Method: hold card at EasyDirect (see EasyDirect Manual)



Technical Data: Card type NXP ICODE SLIX, Passiv RFID HF 13.56MHz, size 1024 bit, according to ISO15693.